

# INSTRUCTION MANUAL



## **LLG-uniSPEC 2**

Please read the User Manual carefully before use, and follow all operating and safety instructions!

<b>1. Preface</b>	
1. Safety . . . . .	3
2. General . . . . .	3
3. Electrical . . . . .	4
4. Warning . . . . .	4
5. Performance . . . . .	4
6. Radio Interference . . . . .	4
<b>2. Introduction of the Instrument</b>	
1. Brief Introduction of the Instrument. . . . .	5
2. Working Principle . . . . .	6
3. Structure. . . . .	7
4. Unpacking Instructions . . . . .	7
5. Specifications. . . . .	8
<b>3. Install Instrument</b>	
1. Environment Required. . . . .	9
2. Voltage Confirmation . . . . .	9
3. Installation . . . . .	10
<b>4 Operation Introduction</b>	
1. Software Structure. . . . .	11
1.1. Menu Tree . . . . .	11
1.2. Firmware Functions . . . . .	12
1.2.1 Basic Mode . . . . .	12
1.2.2 Quantitative. . . . .	12
1.2.3 Kinetics. . . . .	12
1.2.4 Setup . . . . .	12
2. Basic Operation . . . . .	12
2.1 Choose Photometric Mode . . . . .	12
2.2 Set Wavelength . . . . .	12
2.3 Set Parameters . . . . .	12
2.4 Set the position of auto-cell holder. . . . .	12
2.5 Delete the Input Value . . . . .	12
2.6 Delete the test result and stored data . . . . .	12
2.7 Calibrate 100%T/0Abs. . . . .	12
2.8 Measure Samples . . . . .	12
2.9 Print the test result . . . . .	12
3. Preparation before test . . . . .	13
3.1 Switch on the instrument and begin the system self-test . . . . .	13
3.2 Pre-warm . . . . .	13
3.3 Check the cuvettes. . . . .	13
4. Measurement. . . . .	13
4.1 Basic Mode . . . . .	13

4.2	Quantitative. . . . .	15
4.3	Kinetics. . . . .	21
4.4	Setup . . . . .	24
	4.4.1 Power On/Off D2 Lamp. . . . .	24
	4.4.2 Power On/Off W Lamp . . . . .	25
	4.4.3 Set Date and time . . . . .	25
	4.4.4 Calibrate Dark Current . . . . .	26
	4.4.5 Wavelength Calibration . . . . .	26
	4.4.6 Input the lamps changing Wavelength Point . .	27
	4.4.7 Reload Default. . . . .	26
	4.4.8 Version. . . . .	28

## 5 Instrument Maintenance

1.	Daily Maintenance . . . . .	28
2.	Trouble shooting . . . . .	29
3.	Spare part Replacement . . . . .	30

## 1. Preface

### 1. Safety

The safety statements in this manual comply with the requirements of the HEALTH AND SAFETY AT WORK ACT, 1974.

Read the following before installing and using the instrument and its accessories. This instrument should be operated by appropriate laboratory technicians.

### 2. General

The apparatus described in this manual is designed to be used by properly trained personnel in a suitable equipped laboratory. For the correct and safe use of this apparatus it is essential that laboratory personnel follow generally accepted safe procedures in addition to the safety precautions called for in this manual.

The covers on this instrument may be removed for servicing. However, the inside of the power supply unit is a hazardous area and its cover should not be removed under

any circumstances. There are no serviceable components inside this power supply unit. For this instrument, avoid touching the high voltage power supply at all times.

Some of the chemicals used in spectrophotometry are corrosive and/ or inflammable and samples may be radioactive, toxic, or potentially infective. Care should be taken to follow the normal laboratory procedures for handling chemicals and samples.

### **3. Electrical**

Before switching on the apparatus, make sure it is set to the voltage of the local power supply (**see Installation**).

The power cord shall be inserted in a socket provided with a protective earth contact.

The protective action must not be negated by the use of an extension cord without a protective conductor.

### **4. Warning**

Any interruption of the protective conductor inside or outside the apparatus or disconnection of the protective earth terminal is likely to make the apparatus dangerous. Intentional interruption is prohibited.

Whenever it is likely that the protection has been impaired, the apparatus shall be made inoperative and be secured against any unintended operation.

#### **Note**

NEVER touch or handle the power supply on this instrument due to the high voltage!

The protection is likely to be impaired if, for example, the apparatus

- Shows visible damage
- Fails to perform the intended measurements
- Has been subjected to prolonged storage under unfavorable conditions
- Has been subjected to severe transport stresses

### **5. Performance**

To ensure that the instrument is working within its specification, especially when making measurements of an important nature, carry out performance checks with

particular reference to wavelength and absorbance accuracy. Performance checks are detailed in this manual.

### **6. Radio Interference**

For compliance with the EMC standards referred to in the EC Declaration of Conformity, it is necessary that only shielded

cables supplied by us are used when connecting the instrument to computers and accessories.

## 2. Introduction of the Instrument

### 1. Brief Introduction of the Instrument

1. Brief Introduction of the Instrument This instrument is a single beam, general purpose instrument designed to meet the needs of the Conventional Laboratory, This instrument is ideal for various applications, such as: Chemistry, Biochemistry, Petrochemistry, Environmental Protection, Food and Beverage Labs, Water and Waste Water Labs and other fields of quality control and research.

This instrument incorporates a 128 x 64 dots matrix LCD display for photometric results, easy operation and wavelength range of 190 nm to 1100 nm. This instrument is ideal for measurements in the visible and ultraviolet wavelength region of the electromagnetic spectrum.

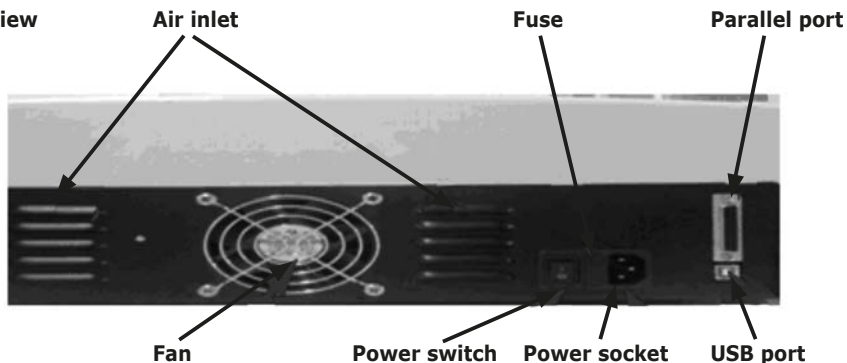


Fig Front View



Fig Control Panel

Fig Back View



## 2. Working Principle

Different matter has different but special absorbance wavelength point. Also, when at the fixed wavelength point, the absorbance has some relation to the substance's (Always transparent Solution) concentration and its thickness. The relation can be concluded as the following Formula which is called Lambert-Beer Law.

$$T = I/I_0$$

$$A = KCL = -\log I/I_0$$

A Absorbance

C Concentration of the Solution

K Absorbance Coefficient of the Solution

L The length of the Solution in the light path

I The intensity of the light focused on the A/D after it permeates the solution to be measured.

$I_0$  The intensity of the light focused on the A/D before it permeates the solution.

### **Note:**


When test, the solvent is usually taken as the Reference Solution and its Transmittance is considered as 100 %T. While the Transmittance of the sample to be tested is a relative value which is got comparing to that of the Reference.

### 3. Structure

The spectrophotometer consists of five parts:

1. Halogen or deuterium lamps to supply the light
2. A Monochromator to isolate the wavelength of interest and eliminate the unwanted second order radiation
3. A sample compartment to accommodate the sample solution
4. Detector to receive the transmitted light and convert it to an electrical signal; and A digital display to indicate absorbance or transmittance. The block diagram below illustrates the relationship between these parts.

Fig **Block diagram for the Spectrophotometer**

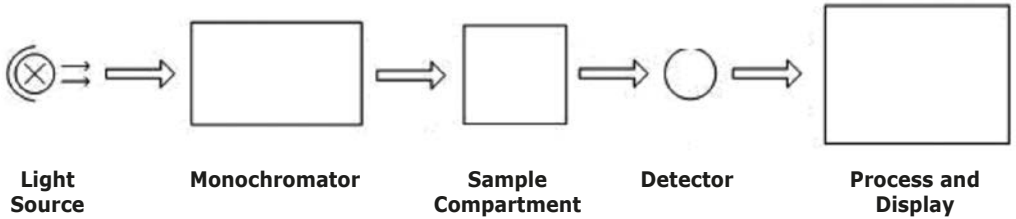

In your spectrophotometer, light from the lamp is focused on the entrance slit of the monochromator where the collimating mirror directs the beam onto the grating. The grating disperses the light beam to produce the spectrum, a portion of which is focused on the exit slit of the monochromator by a collimating mirror. From here the beam is

passed to a sample compartment through one of the filters, which helps to eliminate unwanted second order radiation from the diffraction grating. Upon leaving the sample compartment, the beam is passed to the silicon photodiode detector and causes the detector to produce an electrical signal that is displayed on the digital display.

### 4. Unpacking Instructions

Carefully unpack the contents and check the materials against the following packing list to ensure that you have received everything in good condition.

<b>Packing List Description</b>	<b>Quantity</b>
Spectrophotometer .....	1
Power Cord .....	1
Glass Cuvettes .....	1 Set of 4
Quartz Cuvettes .....	1 Set of 2
Operation Manual .....	1

**Note:**

The Printer and Auto-cell Holder mentioned in this manual are all optional accessories, they do not come standard with the instrument.

## 5. Specifications

Model	LLG-uniSpec 2
Wavelength Range	190-1100 nm
Band Width	2 nm
Wavelength Accuracy	± 0.5 nm
Wavelength Repeatability	0.3 nm
Wavelength Setting	Auto
Photometric Accuracy	±0.5 % T
Photometric Repeatability	0.3 % T
Photometric Display Range	0-200 % T, -0.3-3.0 A
Stability	0.002 A/h @ 500 nm
Stray Light	≤0.05 % T @ 220 nm, 360 nm
Data Output Port	USB
Printer Port	Parallel Port
Display	128 x 64 Dots LCD
Lamps	Deuterium Lamp & Tungsten Halogen Lamp
Detector	Silicon Photodiode
Power Requirement	220 V/50 Hz or 110 V/60 Hz
Dimension (L x W x H)	460 x 380 x 180mm
Net Weight	18 kg



## 3. Install Instrument

### 1. Environment Required

To ensure the best performance, the following conditions are required:

- The best working temperature range is 16-30 °C and the humidity is 45-80 %.
- Keep it as far as possible away from the strong magnetic or electrical fields or any electrical device that may generate high-frequency fields.
- Set the unit up in an area that is free of dust, corrosive gases and strong vibrations.
- Remove any obstructions or materials that could hinder the flow of air under and around the instrument.
- The power requirement is 220±22V@50±1Hz or 110±11V@60±1 Hz.
- Use the appropriate power cord and plug into a grounded outlet.
- If the local voltage is not stable enough, a manostat is required.
- Be away from direct sunlight

### 2. Voltage Confirmation




**Be sure to set the instrument's Voltage Switch at your local power supply, or severe damage may occur! The Voltage Switch is on the bottom of the instrument.**

e.g.: If your local Voltage is 220V, you should set the Voltage Switch at the following status.



If your local Voltage is 110V, you should stir the switch to the Left way by finger, so that it can be set at 110V.

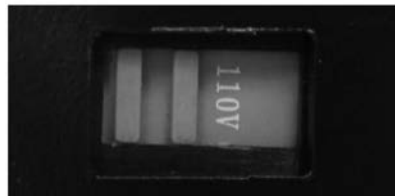

### **3. Installation**

#### **Step 1: Check the packing list**

Unpack the contents, check the materials with the Packing List. Any damage or Lost found, please contact us or the local dealer.

#### **Step 2: Position**

Place the instrument on the stable table carefully.

#### **Step 3: Install printer (Optional)**

Make sure the printer is power off; link the printer's data cable to the instrument's parallel port.

#### **Step 4: Link the power cord**

Make sure the instrument's power switch is in the Off condition, link the power cord to the instrument and insert another end in the socket provided with a protective earth contact.


#### **Step 5: Switch on the power**

Check again, make sure that all the links are right. Switch on the power. The instrument can be used to measure after self-test and pre-warm.

## 4. Operation Introduction

### 1. Software Structure

#### 1.1 Menu Tree

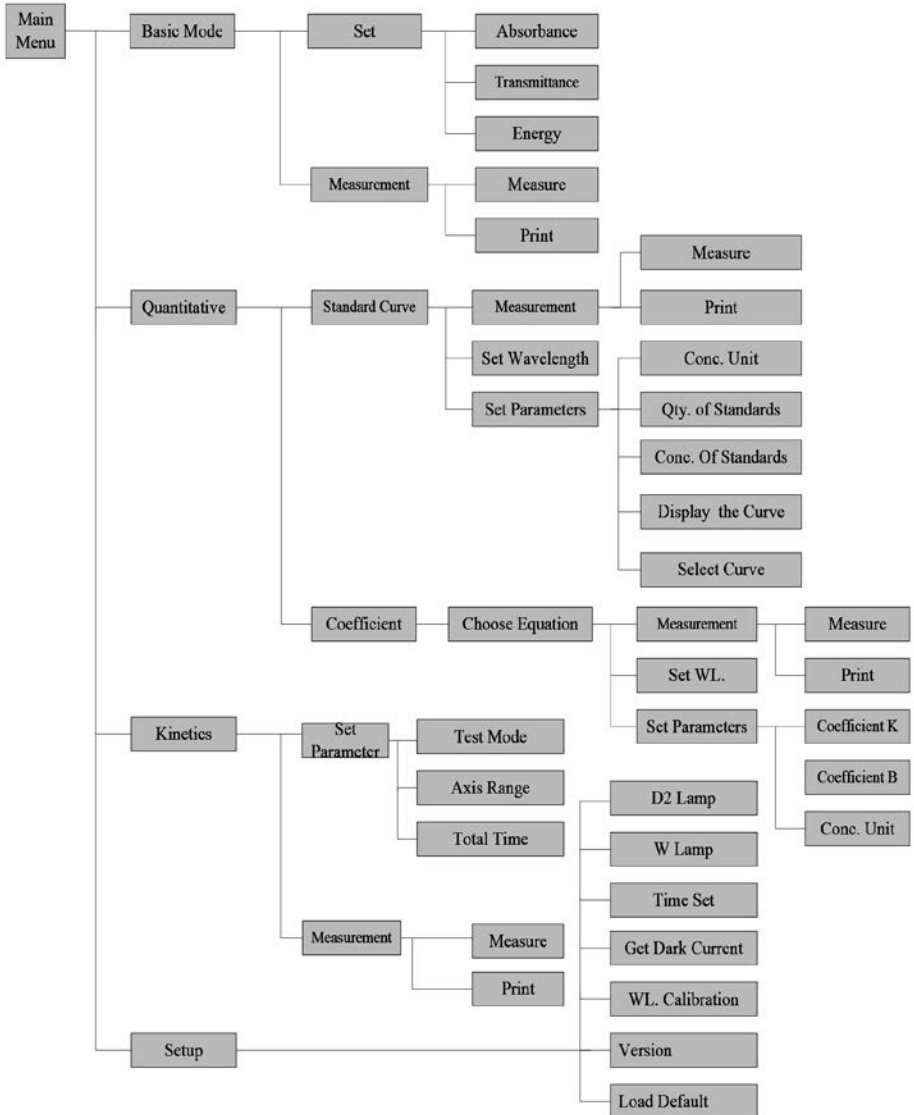

## **1.2 Firmware Functions**

The firmware consists of 4 functions: Basic Mode, Quantitative, Kinetics and System Functions.

### **1.2.1 Basic Mode**

Measure the Absorbance, Transmittance or Energy in a fixed wavelength point. The test result can be stored in the RAM.

### **1.2.2 Quantitative**

**Standard Curve Method:** Set up standard curve by standard samples; use the new curve to measure the unknown sample's concentration. The curves and the test results can be stored in the RAM.

**Coefficient Method:** Input the values of the curve equation's coefficient, and then measure the unknown sample's concentration.

### **1.2.3 Kinetics**

Measure the absorbance or transmittance of a sample at a specific wavelength in a certain time range, the curve can display on the screen and the data can be printed out.

### **1.2.4 Setup**

Light source management; Get Dark Current; Calibrate Wavelength and Set time, etc.

## **2. Basic Operation**

### **2.1 Choose Photometric Mode**

Press Arrow key to choose and  to confirm.

### **2.2 Set Wavelength**

In the measurement Interface, press  to set wavelength, use numeric keyboard to input the value, press  to confirm and set 100%T/0Abs automatically.

### **2.3 Set Parameters**

Press  to go into the parameter setting interface, press Arrow Key to choose the corresponding menu and use numeric key to input the parameter, press  to confirm, press  to return.

### **2.4 Set the position of auto-cell holder.**

(The Auto Cell Holder is optional)

In the measurement Interface, press  first, then press the numeric key of 1-8, and then the corresponding cell will be positioned at the light path.

### **2.5 Delete the Input Value**

Press  to delete the last character of the input.

### **2.6 Delete the test result and stored data**

In the measurement Interface, press  to delete the test result and stored data.

### **2.7 Calibrate 100%T/0Abs**

Put the reference in the light path, press  to calibrate 100%T/0Abs.

### **2.8 Measure the samples**

Put the samples in the light path and press  to measure.

### **2.9 Print the test result**

In the Measurement Interface, press  to print the result

### 3. Preparation before test

#### **3.1 Switch on the instrument and begin the system self-test (Diagnostic Program)**

Make sure to remove all the blocks in the light path; Close the compartment lid and switch on the power; Then the system will begin self-test.

#### **Note:**

During the Diagnostic Program runs, please DON'T open the lid of the compartment!

#### **3.2 Pre-warm**

When the diagnostic program finished, it goes into pre-warm condition. 20 minutes is required before measure.

#### **3.3 Check the Cuvettes**


The cuvette should be clear and there's no solution remains on the transparent surface. If the wavelength is below 340 nm, Silicon Cuvette is a must.

### 4. Measurement

#### **4.1 Basic Mode**


##### **Step 1: Go into Basic Mode interface**

In the main menu, use Arrow Key to choose "Basic Mode", press  to go into corresponding setting interface.

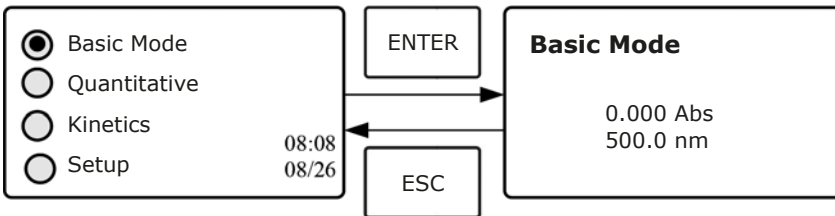


##### **Step 2: Set Wavelength**

Press  to set Wavelength, input the wavelength value by the numeric Keypad, then press  to confirm.

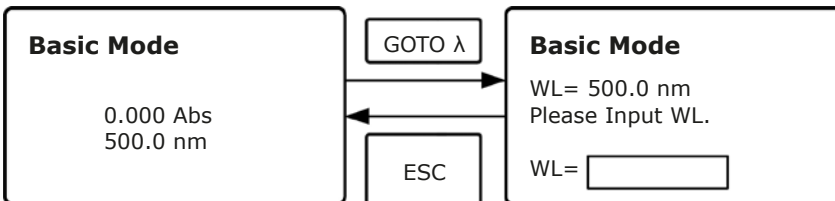


### Step 3: Set Photometric Mode

Press **SET** to go into Parameter Setting Interface, press Array Key to select "Abs", "%T" or "Energy" mode, press **ENTER** to confirm. Press **ESC** to return.

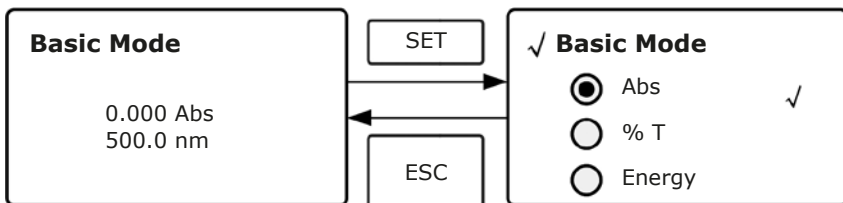

### Step 4: Go into Continuous Measurement Interface

Press **START STOP** to go into continuous measurement interface.

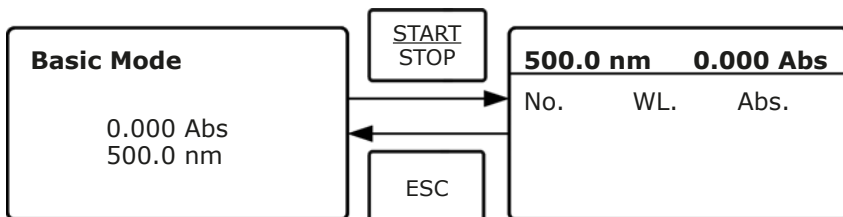

### Step 5: Calibrate 100%T/0 Abs

Put the Reference Solution in the light path, press **ZERO** to calibrate 100 %T/0Abs.

### Step 6: Measure Sample

Put the unknown concentration sample in the light path, press **START STOP** to measure, and the result displays in the data sheet. At the same time, the test result will be stored in the RAM automatically.

Repeat this step to finish all the samples measurement.

### Step 7: Print the data

Press **PRINT** to print the test result.


### Step 8: Delete the data

Press **CLEAR** to delete the test data.

## 4.2 Quantitative


### Step 1: Go into Basic Mode interface

In the main menu, press Arrow Key to choose "Quantitative"; Press  to enter the Quantitative method selecting Interface.

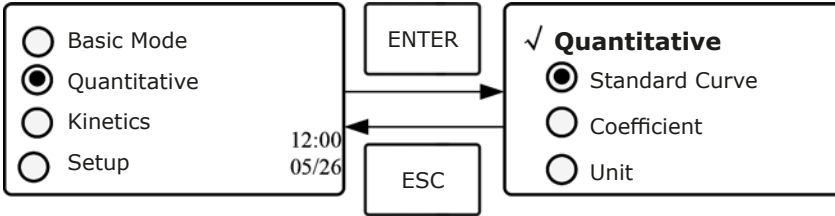

### Step 2: Choose Unit

Move the cursor on "Unit", press  to go into concentration unit choosing interface. Use Arrow Key to move the cursor on the unit you want, and press  to confirm.

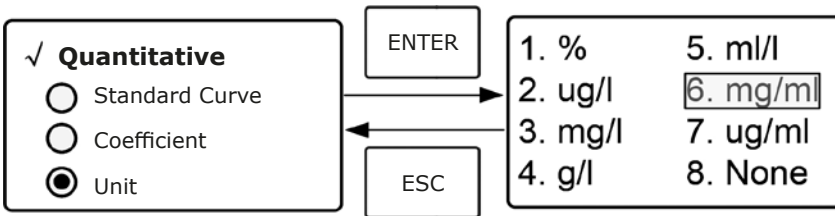


### Step 3: Choose Methods

There are two methods to be selected: Standard Curve and Coefficient.

#### I. Standard Curve:


##### (1) Go into "Standard Curve" Interface

In "Quantitative" menu, press Arrow Key to choose "Standard Curve"; Press  to go into sub menu. Here you can create a new curve, or you can load a saved curve. If you want to delete a saved curve, just choose "Delete Curve".

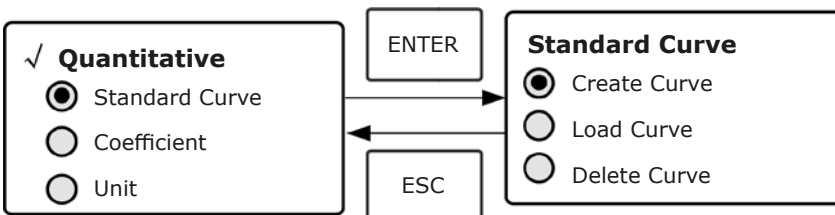


## (2) Set Wavelength


Move the cursor on "Create Curve", then press **ENTER** to go into the Fit Method select interface. If you select "Linear Zero", then the curve must go through the zero point. If you select "Linear Fit", then it doesn't necessarily go through zero point.\*

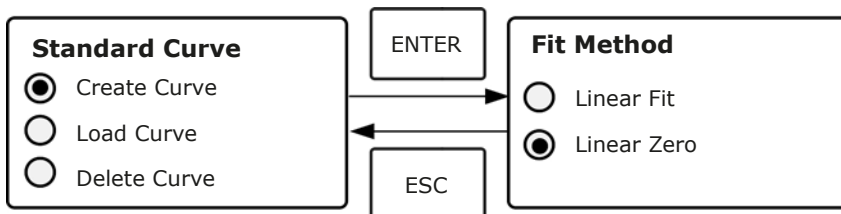

\* After Fit method selected, press **ENTER** to go into the following interface.

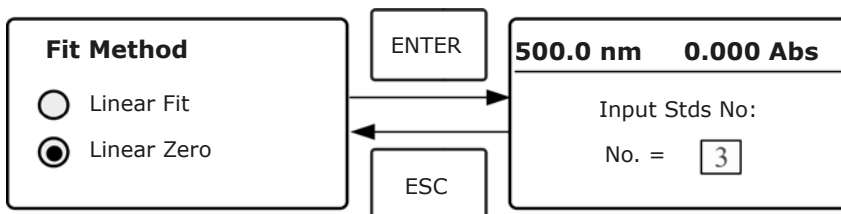

\* Press **GOTO λ** to go into the wavelength setting interface, input the WL. Value by the numeric keypad and press **ENTER** to confirm.

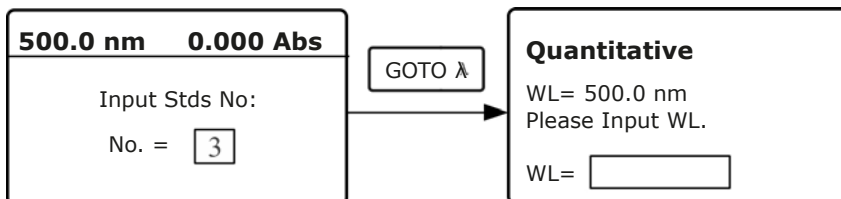


## (3) Calibrate 100 %T/0Abs

Put the Reference Solution in the light path, then press **ZERO** to calibrate 100 %T/0Abs.




**(4) Confirm the quantity of standard samples**


Use the numeric keypad or Arrow Key to input the quantity of the standard samples (e.g. There are 6 Standard Samples), then press  to confirm. The system will go into the following interface.

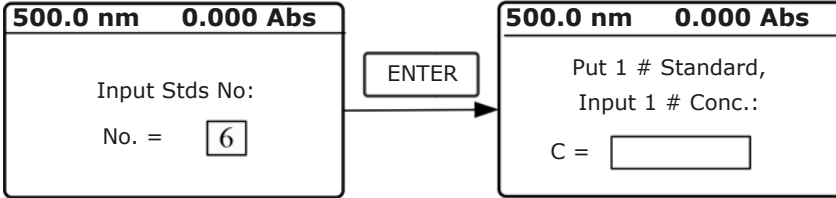**(5) Measure Standard Samples' Concentration**

\* Pull the 1# Standard Sample in the light path and input the value of its concentration by the numeric keypad, then press  to confirm. The system will go into the next interface.

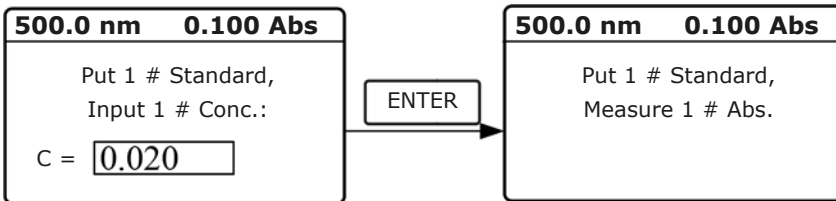


\* Press  to measure 1# Standard Sample, then it goes into the next Interface.

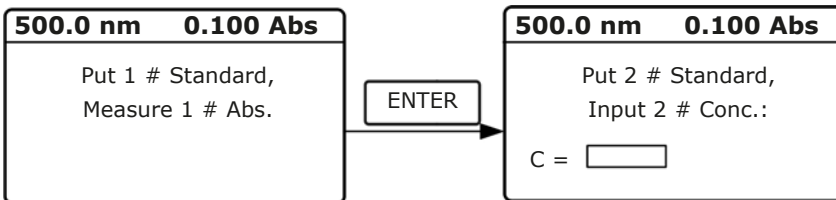

\* Pull the 2# Standard Sample in the light path and input its concentration, then press  to confirm. Press  again to measure 2# sample's concentration.

\* Measure the other standard samples in the same way. When the last sample is finished, press **ENTER** to confirm. Then the Standard Curve and its equation will display on the screen automatically. At the same time, the Curve Equation will be saved in the RAM automatically.

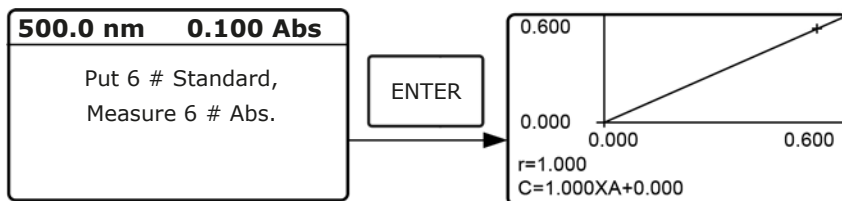


**Note:**

If there's some mistake during the operation, the system will beep 3 times and return the initial interface automatically. Of course, the curve can not be displayed on the screen.

**(6) Measure unknown Samples**

\* Pull the Reference solution in the light path and press **ENTER** to go into the continuous measurement interface.

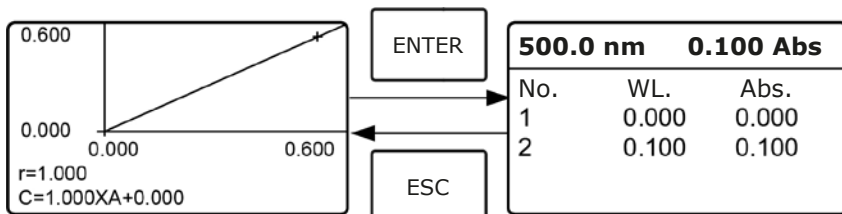


\* Pull the concentration unknown sample in the light path, then press **START STOP** to measure. The result will be displayed on the screen one by one. At the same time, the data will be saved in the RAM automatically. Total 200 groups of data can be saved.


**(7) Load Curve**

All the curve equations will be saved in the RAM automatically. If you want to reload the saved curve, just follow the following steps:

\* Move the cursor on "Load Curve" and press **ENTER** to confirm. Then the system will go into the curve equation record interface.

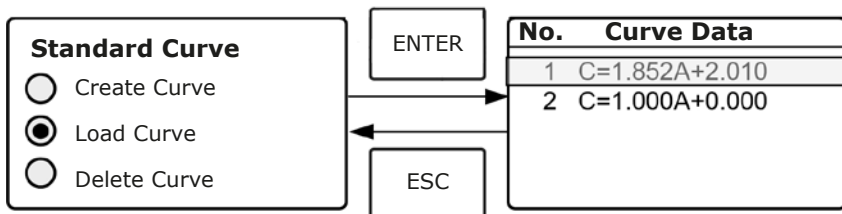

\* Move the cursor on the equation that you need and press **ENTER** to confirm. Then the corresponding curve will display on the screen.

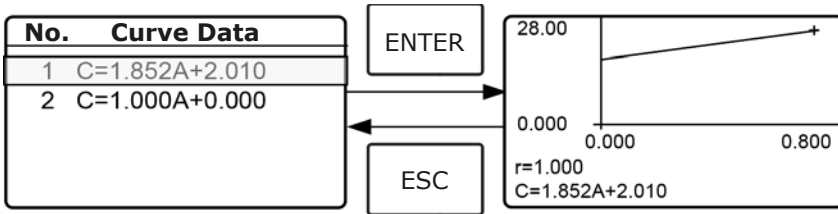


\* Press **ENTER** to go into sample continuous measure interface and measure samples.

**Note:**


When you reload the curves, the wavelength will be set at the point where you create the Standard Curve automatically.

**(8) Delete Curve**

\* Move the cursor on "Delete Curve" and press **ENTER** to confirm. Then the system goes into the curve equation choosing interface.

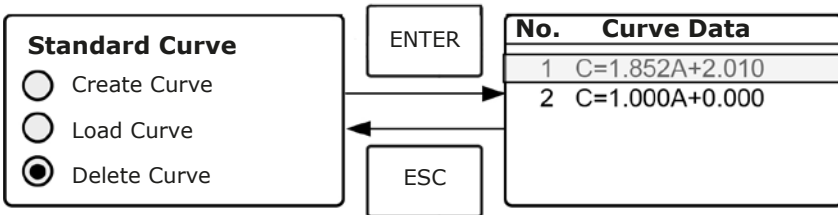

\* Move the cursor on the curve equation that you don't need again and press **ENTER**, then system will ask you to confirm again.

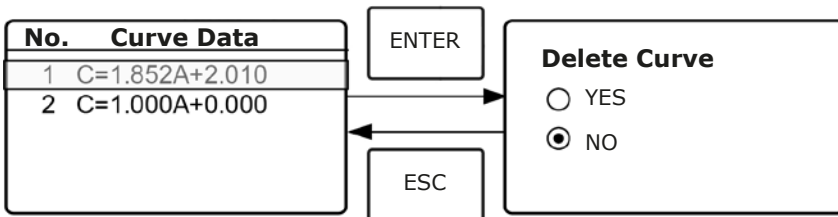


\* Move the cursor on "Yes" and press **ENTER** to confirm, then the curve equation will be deleted. If you don't want to delete it, just choose "NO" or press **ESC** to return.

## II. Coefficient Method:


### (1) Go into "Coefficient" method.

Press Arrow Key to choose "Coefficient", press  to go into pre-test interface, then press  to go into the parameter setting interface.

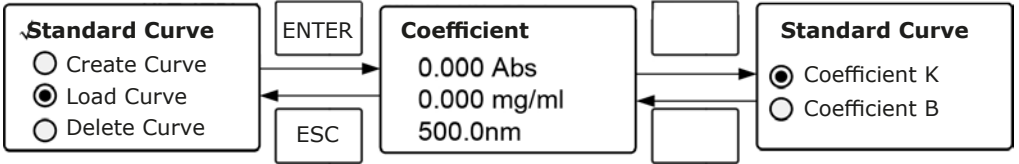

### (2) Parameter Setting


\* Use Arrow Key to move the cursor on "Coefficient K" and press  to go into K setting interface. Input K and press  to confirm, the system will return the former interface automatically.

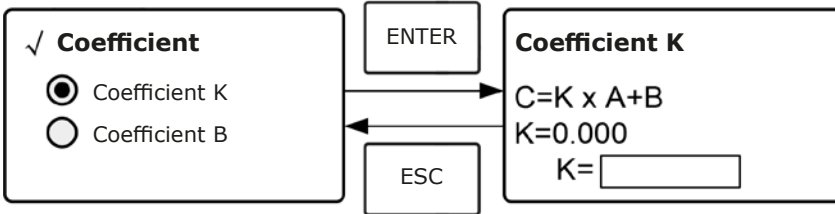

\* Input the value of B in the same way, press  to go back to the pre-test interface.


### (3) Set Wavelength

Press  to go into wavelength setting interface, input the wavelength value by the numeric keypad and press  to confirm.

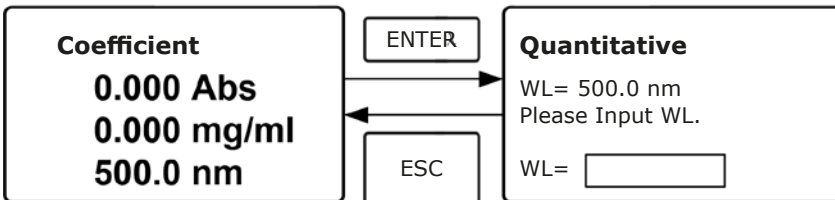

**(4) Calibrate 100 %T/0Abs**

Press **START/STOP** to go into continuous measurement interface. Pull the Reference Solution in the light path, then press **ZERO** to calibrate 100 %T/0Abs.

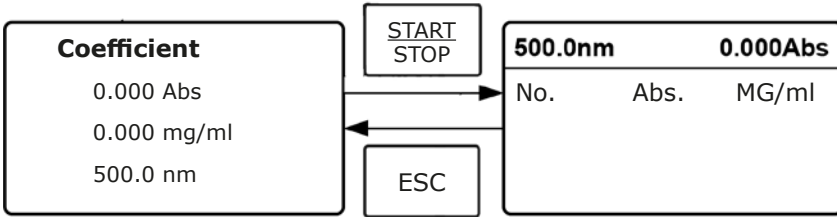

**(5) Measurement**

Place the samples to be measured in the light path, press **START/STOP** to measure, then the result will display on the screen. Repeat this step to measure other samples. All the test results will be listed on the screen and stored in the RAM automatically. They can be reloaded even after a sudden power cut.

**Step 4: Print the test result**

In the continuous measurement interface, press to print the test result.


**Step 5: Delete the test data**

In the continuous measurement interface, press to delete the test result.

**4.3 Kinetics**


This function may be used for time course scanning. The followings are the operation steps for kinetic analysis:

**Step 1: Enter Kinetics measurement interface.**

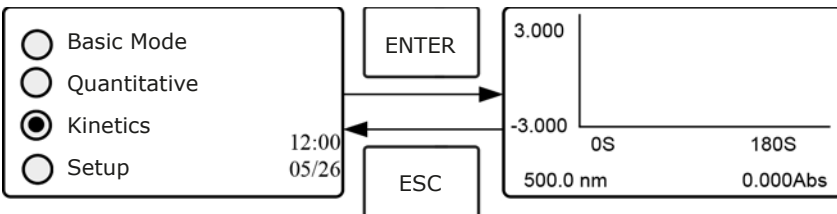


## Step 2: Set parameters

Press  to enter Kinetics setting interface, you can set test mode, kinetics time (Total Time), and the display range of Y-axis (Top Limit and Bottom). You can also browse the test data after completing the measurement by selecting "Record List".

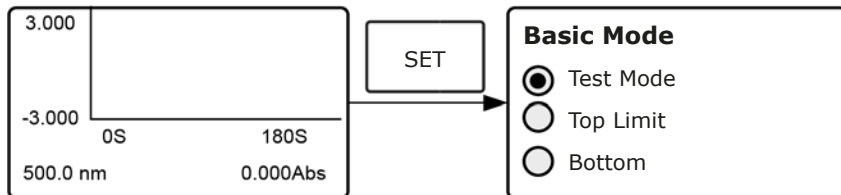


### \* Test mode

There are three test modes for chosen, Absorbance, Transmittance and Energy. Press Arrow Key to select the wanted mode and press  to confirm.

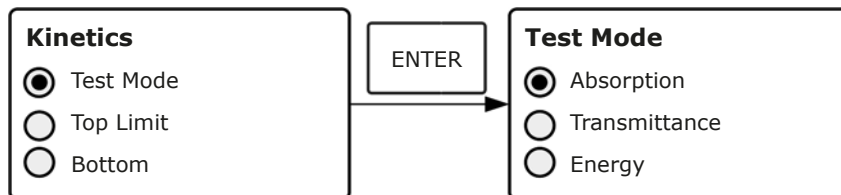


### \* Y-axis Setting

Press Arrow Key to select "Top Limit" and press  to confirm. Input the value according to certain test mode, press  to confirm. You can set the Bottom value in the same way.

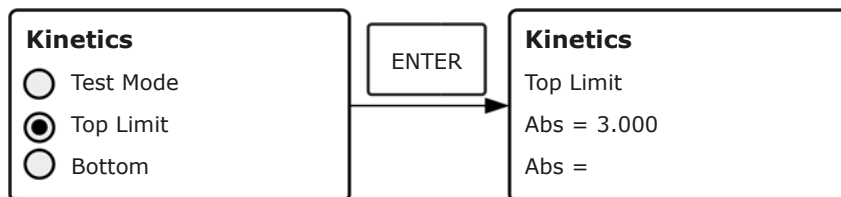

### \* Test time Setting

Press Arrow Key to select "Total Time" in Kinetics setting interface and press  to go into the setting interface. Input the time and press  to confirm.

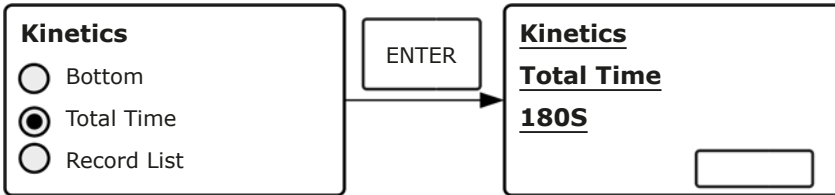


#### Note:

The valid range of total time is 10 ~ 3600.


### Step 3: Set the measurement wavelength.

In Kinetics measurement interface, press  to enter the wavelength setting interface (Fig. 4-30). Input the wavelength value and press  to confirm. When get to the setting wavelength point, it will return to Kinetics measurement interface.

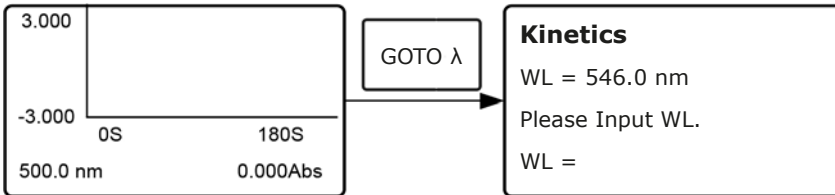

### Step 4: Kinetics measurement

In Kinetics measurement interface, put the blank solution into the light path, press  to adjust blank. Then pull the sample solution in the light path and press , it will start the kinetics measurement and a dynamic curve shows in real-time.

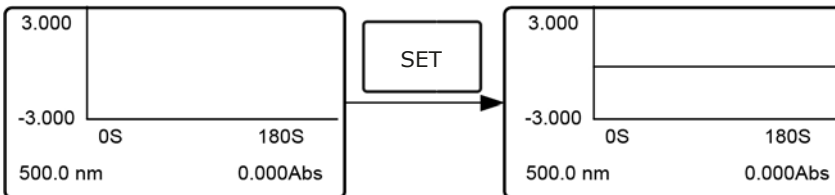


#### Note:

If users want to stop kinetics measurement during the course, just press the key of .


### Step 5 Check the Test Data

When the kinetics measurement finished, you can select Record List to check all the data.

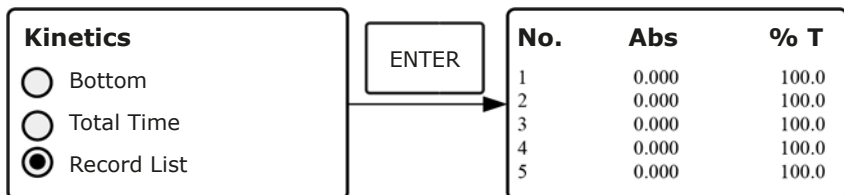

### 4.4 SET UP


In the main menu, press Arrow Key to choose "Setup" and press  to go into utility setting interface.

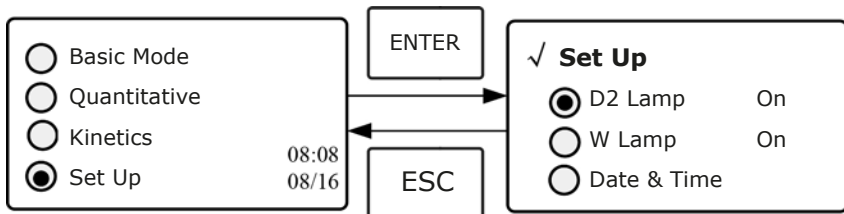

#### 4.4.1 Power on/off D2 lamp

**i** When the wavelength point needed is in the range of 340-1100nm, the D2 lamp can be switched off to prolong its life time.

Use Arrow key to choose "D2 Lamp" and press  to go into D2 lamp setting interface. Press Arrow key to choose "On" or "Off" and press  to confirm. Press  to return.

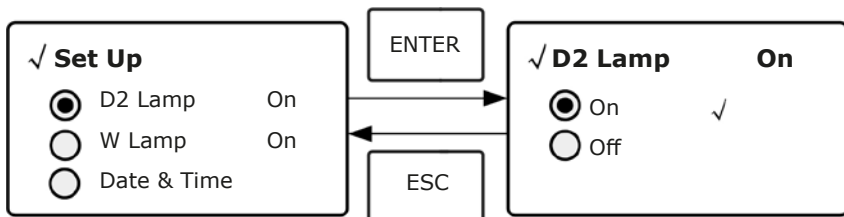




#### 4.4.2 Power on/off W Lamp


**i** When the wavelength point needed is in the range of 190-339 nm, the W lamp can be switched off to prolong its life time.

Use the key of Arrow Key to choose "W Lamp", press  to go into W lamp setting interface. Use the key of Arrow Key to choose "On" or "Off" and press  to confirm. Press  to return.

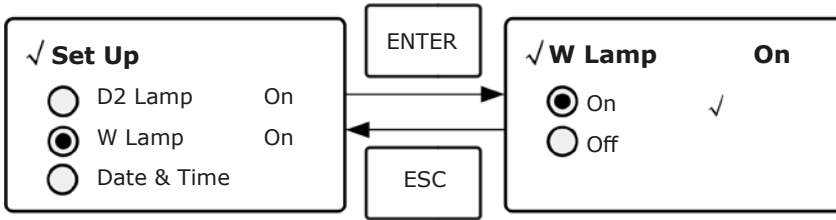

#### 4.4.3 Set Date & Time

Press Arrow Key to choose "Date & Time" and press  to go into setting interface. Use Arrow Key to choose setting item and input the value by numeric keypad. (Hour: 0-23, Minute: 0-59, Second: 0-59, Year: 0000-9999, Month: 1-12, Day: 1-31, Week: 1-7), press  to return.

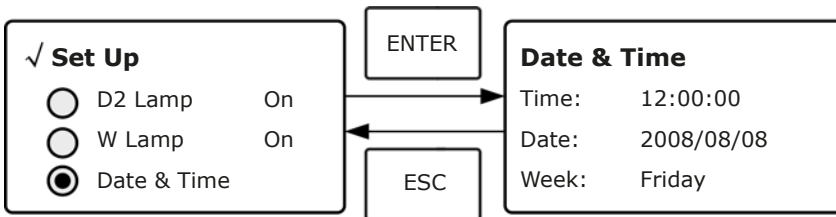

#### 4.4.4 Calibrate Dark Current


**i** When the working condition changes, A dark current calibration is required before any measurement.

Use Arrow key to choose “Dark Current” and press  to confirm.

**Note:**

Remove the cuvettes from the light path before this action, don’t open the lid of the compartment during the course of calibration.


Press  to return when it finishes.

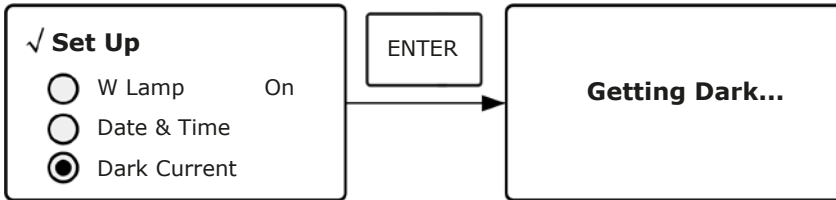

#### 4.4.5 Wavelength Calibration

**i** After some time, the energy of lights would drop off and there will be some but slight influence on the test result. In this case, users can reset the wavelength to compensate. We suggest the users to reset the wavelength every one or two months.

Remove all the blocks in the light path and close the lid of the compartment, press Arrow Key to choose “Calibrate WL.” and press  to begin the calibration (Fig. 4-38). The lid of the compartment should not be opened during the calibration. It returns automatically after calibration.

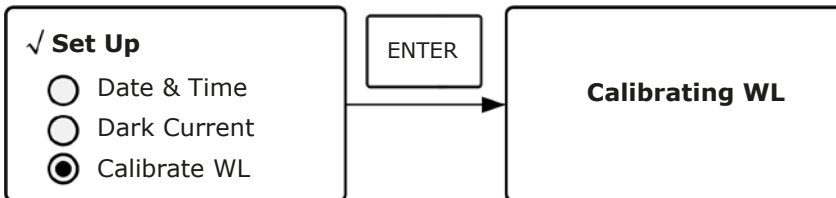


#### 4.4.6 Lamps Change (Input the lamps changing wavelength point)




The instrument permits the users to set the lamps changing wavelength point. In the range of 300-400nm, users can choose the lamp freely.

Press Arrow Key to choose "Lamp Change" and press  to go into corresponding setting interface. Input the value of lamps changing wavelength point by the numeric keypad (300-400 nm), then press  to confirm and return.

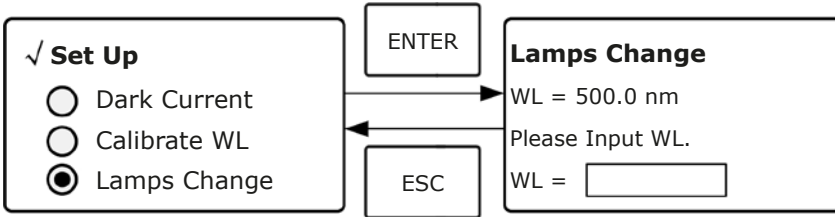


#### 4.4.7 Load Default

Press Arrow Key to choose "Load Default" and press  to confirm. Press Arrow Key to choose "Yes" and press  to confirm. Then the system will recover the default setting when leave factory. Choose "No" with  pressed to cancel the operation and return.

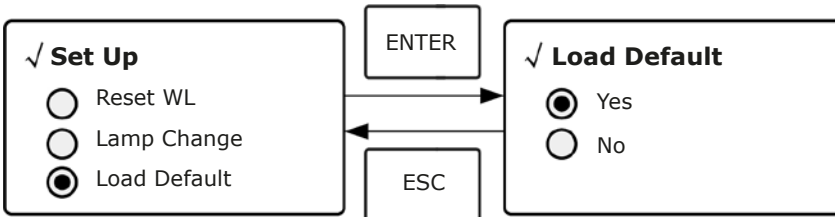

#### 4.4.8 Version

Press Arrow Key to choose "Version" and press  to confirm. You'll see the version of the software and hardware. Press  to return.

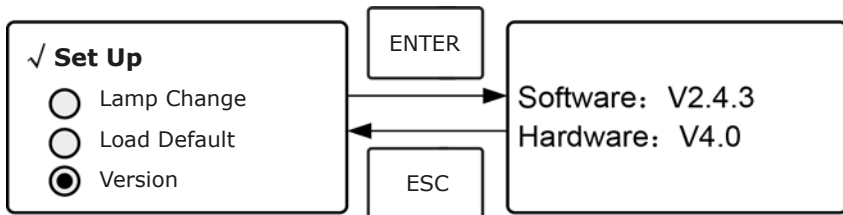

## 5. Instrument Maintenance

To keep the instrument work in good condition, constant maintain is needed.

### 1. Daily Maintain

#### **1.1 Check the compartment**

After measurement, the cuvettes with sample solutions should be taken out of the compartment in time. Or the volatilization of the solution would make the mirror go moldy. Users must pay more attention to the corrosive sample and liquid easy to volatilize. Any solution remains in the compartment should be wiped off immediately.

#### **1.2 Surface Clean**

The cover of the instrument is with paint. Please use wet towel to wipe off the drip on the surface immediately. Organic solution is forbidden to be used to clean the cover. Please wipe off the dirt on the cover timely.

#### **1.3 Clean the cuvettes**

After every test or after a solution change, the cuvettes should be cleaned carefully, or the remains on the surface would cause measuring error.

## 2. Troubleshooting

### 2.1 Dark Current Error when self-test

**Possible Cause**

- Open the lid of the compartment during the course of self-test.

**Solution**

- Close the lid of the compartment and switch on the power again.

### 2.2 No response after power on

**Possible Cause**

- Bad contact in power supply
- Fuse melt

**Solution**

- Improve the contact
- Replace a new fuse

### 2.3 Printer not work, printer error

**Possible Cause**

- No power supply
- Bad contact in power supply
- Bad contact of the data cable

**Solution**

- Switch on the power supply
- Improve the contact
- Improve the contact

### 2.4 No Stable Reading

**Possible Cause**

- No enough pre-warm
- Glass cuvettes used in UV Range
- No stable Sample
- Much higher sample concentration
- Low voltage or unstable power supply
- Lights defect
- Light used up

**Solution**

- Increase the pre-warm time
- Use Silicon Cuvettes
- Improve the sample
- Dilute the sample
- Improve the power condition
- Replace a new lamp
- Replace a new lamp

### 2.5 Worse Repeatability

**Possible Cause**

- Unstable sample
- Cuvettes polluted

**Solution**

- Improve the sample
- Clean the cuvettes

### 2.6 Incorrect reading

**Possible Cause**

- Dark Current Error
- Worse matching of the cuvettes

**Solution**

- Re-get the dark current
- Improve the matching of the cuvettes

### 3. Spare part replacement

#### Replace the Fuse



**Danger! Be sure to switch off the power and unplug the socket before replacement!**

#### Step 1: Tools preparation



Prepare a 3×75 Flat Blade screwdriver

#### Step 2: Switch Off the power supply

Switch off the power supply, and unplug the socket.

#### Step 3: Take out the Fuse Seat

Take out the fuse seat by the screwdriver.

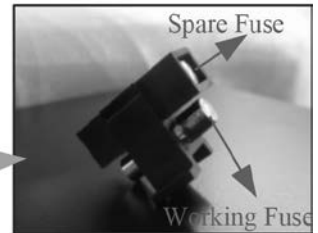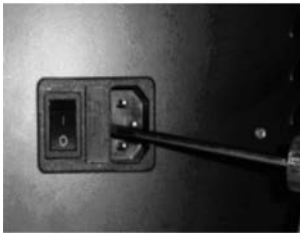

#### Step 4: Replace a new fuse


Pick out the spare fuse and replace it to the working position.

#### Step 5: Reset the fuse seat

Replace the fuse seat in the power socket

#### Step 6: Switch on the power

Plug the socket and switch on the power

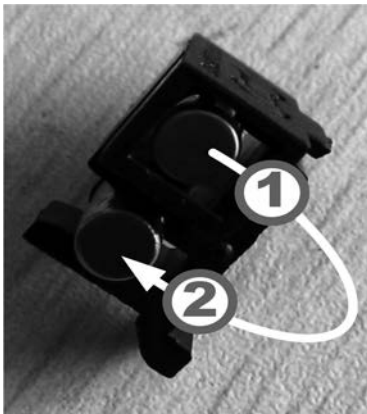

## WARRANTY REPLACEMENT

This product is warranted to be free from defects in material and workmanship for a period of three (3) years from date of purchase.

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in this instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces or other causes not arising from defects in original material or workmanship. This warranty does not cover any incidental or consequential damages, commercial loss or any other damages from the use of this product.

The warranty does not cover damage to paint or finish and defects or damages caused by physical and chemical abuse or normal wear and tear. The warranty is invalidated by any non-factory modification, which will immediately terminate all liabilities on us for the products or damages caused by its use. The buyer and its customer shall be responsible for the product or use of products as well as any supervision required for safety. If requested the products must be returned to the distributor in well packed and insured manner and all shipping charges must be paid.

Information on the disposal of electrical and electronic devices in the European Community:

The disposal of electrical devices is regulated within the European Community by national regulations based on the European Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). According to these regulations, any devices supplied after 13.06.05 in the business to business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. They are marked with the following symbol to indicate this. As disposal regulations within the EU may vary from country to country, please contact your supplier if necessary.

### PRODUCT DISPOSAL



In case the product is to be disposed of, the relevant legal regulations are to be followed.



**Lab Logistics Group GmbH**  
Am Hambuch 1  
D-53340 Meckenheim/Deutschland

Tel.: +49 (0)2225 9211- 0

Fax: +49 (0)2225 9211-11

[www.llg-labware.com](http://www.llg-labware.com)

[info@llg-labware.com](mailto:info@llg-labware.com)