Operations Manual LLG-uniCFUGE 2

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Introduction:

Thank you for purchasing LLG-uniCFUGE 2. Please read this manual thoroughly prior to operating the instrument.

Specifications LLG-uniCFUGE 2:

| | |
|-------------------|------------------------|
| Speed | 4000 rpm & 6000 rpm |
| RCF | 1200 g & 2000 g |
| Dimension (DxWxH) | 135x147x120 mm |
| Power | 100-240 V~ 50-60 Hz 1A |
| Weight | 0.7 kg |

The rotor may only be operated on this unit with the specified speeds. The device is not suitable for operation in humid environments and must not be autoclaved.

Switchable Speed options (6K or 4K)

Set the speed switch on the bottom of the device, then select the speed you need, 6K for 6000 rpm and 4K for 4000 rpm.

LLG-uniCFUGE 2 is supplied with 1 rotor:

- 6 x 1,5/2,0ml , 6 x 0,5ml tubes and 2-8 x 0,2ml PCR Strip rotor

Rotor Exchange for LLG-uniCFUGE 2:

- **Removing the Rotor from the motor shaft:**
Use the included allen key to loose the set screw of the rotor by turning counter-clockwise 4-5 complete turns. The rotor can now be removed by lifting it evenly off of the rotor shaft.
- **Attaching the Rotor to the motor shaft:**
Place the rotor on the motor shaft and press down until the rotor comes to a stop. The set screw can now be tightened by inserting the allen key and turning clockwise 4-5 complete turns.

Operation:

- Attach the power cord and place the power switch into the on "-" position.
- Open the lid and load your samples into the rotor. Always ensure a balanced load (Please see the following section, **Loading the Rotor**, for details.)
- Close the lid, the rotor quickly accelerates to the maximum speed.
- When the desired run time is achieved, push on the lid button to open the lid, the rotor stops and the samples can be retrieved.

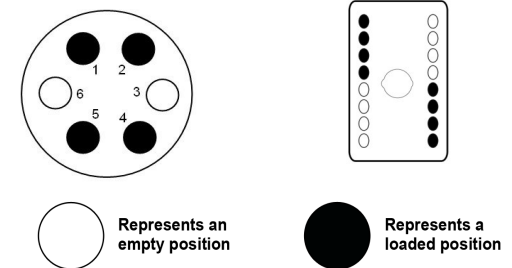
NEVER Attempt to remove samples before the rotor has come to a complete stop.


Loading the Rotor:

- To ensure safe operation and long life of the device, samples must always be loaded into the rotor in a balanced manner.
- Tubes and strips must always be loaded symmetrically to each other and contain the same volume of liquid. See below:

Examples of properly balanced rotors:

Examples of improperly balanced rotors:

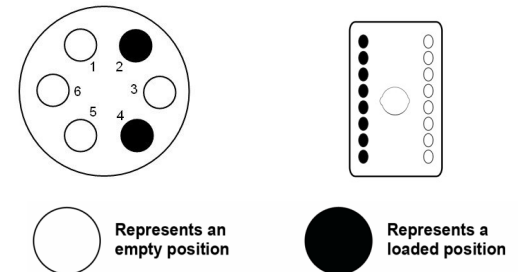Cleaning and Maintenance:

The centrifuge rotors can be removed for cleaning by using the included allen wrench to loosen the set screw. Once removed, the rotors can be cleaned with isopropyl alcohol or can be sterilized in an autoclave at 121°C for 20 minutes.

**Lab Logistics Group GmbH / Am Hambuch 1
D-53340 Meckenheim / www.llg-labware.com**

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