INSTRUCTION MANUAL	Page 4
GEBRAUCHSANWEISUNG	Page 11
MANUEL D'INSTRUCTIONS	Page 17
MANUAL DE INSTRUCCIONES	Page 23

## **LLG-uniCFUGE 3 pro**

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Thank you very much for purchasing this mini Centrifuge LLG-uniCFUGE 3 pro.

Please read the following instructions carefully and always keep this manual within easy reach.

<b>1. Introduction</b>	<b>4</b>
<b>2. Intended Use</b>	<b>4</b>
<b>3. Features</b>	<b>4</b>
<b>4. Standard Accessories</b>	<b>4</b>
<b>5. Technical Specifications</b>	<b>5</b>
<b>6. Safety Precautions</b>	<b>5</b>
<b>7. Installation</b>	<b>6</b>
<b>8. Rotor Balancing</b>	<b>6-7</b>
<b>9. Interchangeable Rotor</b>	<b>7</b>
<b>10. Changing the Rotor</b>	<b>7</b>
<b>11. Operation</b>	<b>7</b>
<b>12. Warranty</b>	<b>8</b>

## Ordering Information

<b>Cat. No.</b>	<b>Description</b>
6.263 522	LLG-uniCFUGE 3 pro, Centrifuge with exchangeable rotor and EU plug
6.263 523	Rotor for 8 x 1,5/2 ml tubes for LLG-uniCFUGE 3 pro
6.263 524	Rotor for 16 x 0,2 ml tubes for LLG-uniCFUGE 3 pro
6.263 527	LLG-uniCFUGE 3 pro, Centrifuge with exchangeable rotor and UK plug

## 1. Introduction

This mini centrifuge offers versatility, simplicity, safety, and ease of usage. The mini centrifuge LLG-uniCFUGE 3 pro is especially suitable for microfiltration, cell separations, centrifugal tube sedimentation, and spin down applications. Please read this manual thoroughly before operating the centrifuge.

## 2. Intended Use

This is used in various laboratories to separate particles from suspension or even macromolecules from solutions according to their density. The different biological substances that are usually separated by centrifugation are microbial cells, organelles, proteins, DNA, and RNA.

**NOTE:** *This user guide is intended to assist with the operation and care of the unit only and not for its repair. The customer should not attempt to service or repair the unit. If repairs are required, please contact your local LLG-Labware distributor.*

## 3. Features

- Mini centrifuge for 8 tubes
- Digital display with micro processor control
- Set variable speed & timer
- Electronic safety brake on lid opening for user safety
- Magnetic rotor for easy tool free changing of rotor
- Starts and stops in seconds
- Lid safety interlock
- Smooth and silent, discreet running
- Small footprint
- Interchangeable magnetic rotors (MCT rotor & PCR rotor)
- Low voltage power consuming device

## 4. Standard Accessories

- 1.5/2 ml rotor, PCR strip rotor
- Adaptor for 0.5 ml tubes, pack of 8
- Adaptor for 0.1/0.2 ml tubes, pack of 8
- Power Adaptor: Inputvoltage: 100 - 240 VAC, 50/60 Hz  
Outputvoltage: 24 V, 1.5 A
- Product User Manual, Warranty Card and Declaration Conformity

## 5. Technical Specifications


Motor Type	BLDC Motor
Rotor Capacity	8 x 1.5 /2.0 ml Tube Rotor & 2 Place PCR Strip Rotor
Speed	1000 - 6000 RPM
	2000 x g (MAX RCF)
Speed Accuracy	± 100 RPM
Time	1 - 25 min. or Infinite
Size (W X D X H)	162 x 157 x 116 mm
Weight	1.1 kg
Input Voltage	24 V, 0.62 A
Power Consumption	15 W

## 6. Safety Precautions

The personal mini centrifuge is designed to be safe under the following conditions when operated with necessary precautions mentioned below:

- Indoor use only.
- Ambient temperature of 4 °C to 45 °C.
- Main supply voltage fluctuations not to exceed +/- 10% of the nominal voltage.
- Position at least 20 cm from walls, 20 cm from ceiling and 20 cm between other products.
- Make sure the rotor is placed in the shaft groove properly. If the rotor is not installed properly, the machine will make noise, turn off the machine & install the rotor as instructed for fitment of the rotor.
- Always load the rotor symmetrically. Each tube should be counterbalanced by another tube.
- Make sure to always use only specified mini centrifuge tubes made from plastic and designed to withstand centrifugal forces of at least 2000 g.
- Do not operate the centrifuge without the rotor being properly attached to the shaft. Re-check by trying to pull it up manually, especially after the rotor has been changed or re-assembled.
- Do not move the centrifuge while it's under operation.
- Do not fill the tubes while they are in the rotor. Liquid spillage may harm the unit.
- Do not put hands in the rotor chamber unless the rotor has completely stopped.
- Do not centrifuge flammable, explosive or corrosive materials.
- Do not use solvents or flammable media near this or other electrical equipment.
- In case the unit is dropped from the table, please check the casing for any cracks/damage.
- The centrifuge and its components are non autoclavable.
- Do not use the centrifuge in any manner not specified in these instructions.
- Use mild soap solution to clean & disinfect. Do not use sanitizer / IPA solutions.

## 7. Installation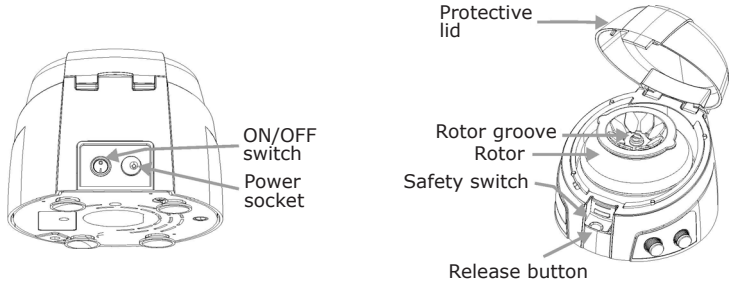

This mini centrifuge is supplied in a packaging box. Open the box, then gently remove the centrifuge from the box and plug in the power adaptor pin before connecting to the main power source. The instruction manual and the accessories should be kept with the centrifuge. Please keep all packaging in safe storage for at least 3 years for warranty purposes. Make sure the rotor is tight to avoid any accident. Load the rotor with the micro tubes for centrifugation. Place the ON/OFF switch to „0“ on back side and close the lid to start centrifugation.

For more clarification refer to the RCF chart below:

RCF CHART (FIXED)					
RPM	ROUND ROTOR (MCT ROTOR)				PCR STRIP
	0.2 ml	0.5 ml	1.5 ml	2 ml	
3000	319	424	468	488	467
4500	718	953	1053	1098	1050
6000	1276	1695	1872	1952	1868


## 8. Rotor Balancing

1. Always balance the rotor before centrifugation. The following are symmetrical loadings of centrifuge tubes to the rotor.

## 8. Rotor Balancing

2. Ensure that the samples in the tubes are of equal volume.
3. Symmetrical loading of tubes in the rotor is necessary, otherwise vibration can occur which can cause serious damage to the centrifuge.
4. The following are the incorrect methods of loading tubes in the centrifuge the rotor.

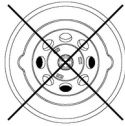## 9. Interchangeable Rotor


Magnetic Rotor with 8 slots  
8 X 0.2, 0.4, 0.5, 1.5, 2 ml

Magnetic PCR Rotor  
16 X 0.2 ml

## 10. Changing the Rotor

This centrifuge features magnetic rotors. Thus changing the rotors can be done without any tool. The rotors fit on the central shaft of the motor through magnet situated at the both the rotors (MCT & PCR rotor). The rotor fits through snap to the central shafts through gentle guidance into the groove on the rotor adapter.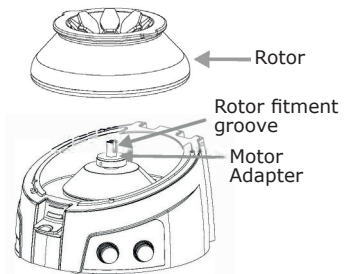

## 11. Operation

To operate the mini centrifuge, the below steps should be followed:

1. Connect the power adaptor & Switch ,On' from the back side.
2. Load the rotor symmetrically to avoid accidents and damage to the centrifuge.
3. Close the lid & flip ON the back side switch to start centrifugation.
4. Rotate\* speed knob to set the speed between 1000-6000 RPM (Display in two digit [x 100] for conversion).
5. Rotate time knob to set time from 1 to 25 min. or infinite.
6. After the rotor has stopped, press the lid release button and open the lid with your thumb in front and fingers on the top - simply lift the lid back on the hinge.
7. If you want to stop the centrifuge quickly, push the lid release button and the rotor will stop within a few seconds.

**\*NOTE:** to increase rotate knob clock wise & to reduce rotate knob counter clockwise.

## 12. Warranty

This product is warranted to be free from defects in material and workmanship for a period of three (3) years from date of purchase.

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in this instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces or other causes not arising from defects in original material or workmanship. This warranty does not cover any incidental or consequential damages, commercial loss or any other damages from the use of this product.

The warranty does not cover damage to paint or finish and defects or damages caused by physical and chemical abuse or normal wear and tear. The warranty is invalidated by any non-factory modification, which will immediately terminate all liabilities on us for the products or damages caused by its use. The buyer and its customer shall be responsible for the product or use of products as well as any supervision required for safety. If requested the products must be returned to the distributor in well packed and insured manner and all shipping charges must be paid.

Information on the disposal of electrical and electronic devices in the European Community:

The disposal of electrical devices is regulated within the European Community by national regulations based on the European Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). According to these regulations, any devices supplied after 13.06.2005 in the business to business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. They are marked with the following symbol to indicate this. As disposal regulations within the EU may vary from country to country, please contact your supplier if necessary.

### PRODUCT DISPOSAL

In case the product is to be disposed of, the relevant legal regulations are to be followed.

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Lab Logistics Group GmbH