## OPERATION MANUAL

# **LLG-uni*INCUBATOR* 20 Cool Cooling Incubator**

Thank you very much for purchasing this  
LLG-uni*INCUBATOR* 20 Cool Cooling Incubator  
Please read the following instructions carefully  
and always keep this manual within easy reach.

## Contents

<b>1. Introduction</b>	<b>3</b>	<b>7. Operation</b>	<b>6</b>
<b>2. Intended Use</b>	<b>3</b>	7.1 Operating the Instrument	<b>6</b>
<b>3. Features</b>	<b>4</b>	7.2 Calibration	<b>7</b>
<b>4. Technical Specifications</b>	<b>4</b>	<b>8. Cleaning</b>	<b>8</b>
<b>5. Installation</b>	<b>5</b>	<b>9. Ordering Information</b>	<b>8</b>
<b>6. Safety Precautions</b>	<b>6</b>	<b>10. Warranty</b>	<b>9</b>

## 1. Introduction

This is a mini cooling incubator that can control temperature in the chamber for incubation experiments. The temperature ranges from -15 °C from ambient to 60 °C. This equipment is of 22 L capacity to hold different flasks up to 2 L.

**Note!**

This device cools based on the ambient temperature of the room. It should not be used outside in the heat. It works best when used indoor in a cool room.

## 2. Intended Use

This instrument is used in order to control temperature and create a homogeneous environment required for cell culture or for biological & non-biological sample incubation. This instrument is intended for indoor use & should only be used by trained laboratory staff.

**Note!**

Before using the instrument, please carefully read this user manual. This user manual is intended to assist with the operation and care of the unit. For repair please contact the supplier.

### 3. Features

- Digital Temperature control from Ambient -15 to 60 °C
- Compact footprint, usable on small bench space
- Accepts multiple types of flask up to 2 L
- Internal power outlet for installing small shaker, mixer or other instruments

### 4. Technical Specifications

<b>LLG oder codes</b>	<b>6.263 572 LLG-uniINCU 20 Cool, EU plug</b> <b>6.263 573 LLG-uniINCU 20 Cool, UK plug</b>
Temperature Range	Ambient -15 °C to 60 °C
Temperature Accuracy	0.5 °C (at 37 °C)
Temperature Uniformity	+/- 1.5 °C (at 37 °C)
Temperature Increments	+/- 1 °C
Exterior Dimension	410 x 340 x 480 mm
Interior Dimension	260 x 235 x 360 mm
Pollution Degree	2
Environment	For indoor use only
Capacity	22 L
Weight	7.4 Kg
Electrical	100-240 V, 50-60 Hz

## 5. Installation

- Place the incubator on a clean, level & stable surface. Position the appliance and allow it to stand for two or three hours (ten hours in summer) before it is turned on for the first time. Do not position the unit next to any heat source e.g. boiler or direct sunlight.
- A clearance of approximately 13-15 cm must be left between the sides of the unit and any adjacent cabinets or walls to allow for adequate air circulation. The unit is not designed for recessed installation.
- Remove the power supply from the white box inside the chamber and plug the wall end into the wall outlet. Plug the other end of the power supply into the 2 pin inlet (top inlet) labelled DC onto the backside of the unit.

### Internal Outlet

To use the internal outlet plug the included power cable into the bottom power inlet on the backside of the instrument, internal outlet is now powered.

**Note!**

The voltage of the internal outlet is dependent on the wall voltage (if the wall voltage is 230 V, voltage of the internal outlet is also 230V).

## 6. Safety Precautions

- Do not use this item near water where it can be exposed to getting wet.
- Make sure the item is only operated from the type of power source indicated on the label.
- Do not add any attachments that were not originally included with the item as they may cause harm to the machine and become a safety issue.
- Make sure the power cords are placed in a position where they will not be damaged.
- Make sure not to overload power outlets by plugging in too many devices in one socket.
- Unplug the unit when not in use for a long period of time.

## 7. Operation


### 7.1 Operating the Instrument

1. Open front of the unit and place samples inside.
2. Plug the adaptor into the machines and then put the electric plug into the wall.
3. Turn power switch to „ON“ or „COOL“.
4. Set the desired temperature ranging from -15 °C from ambient (can go as low up to 2 °C) to 60 °C.
5. The temperature can be changed by pressing „UP“ for increment & „DOWN“ for decrement. The changed values are saved automatically.

## 7.2 Calibration

### Adjustment of the temperature off set

1. Press up and down button simultaneously and hold down until display shows "00A" as the image below.

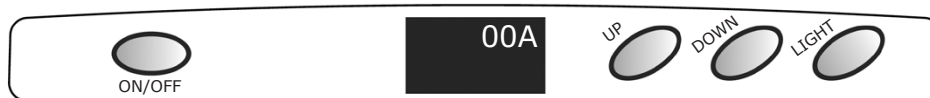2. Set the required value within the range of  $\pm 5$  °C, acknowledge the entry and wait for automatic storage of the value.  
 3. The new temperature is now adjusted.

### For example

- The display value is 30 °C but actual temperature is 28 °C as measured by reference temperature probe.
- Press up and down button simultaneously and hold down together until calibration mode is displayed 00A on the digital display.
- Press down button twice and stop when digital display shows "002A".
- The controller adds the adjusted offset to the displayed temperature.

#### Note!

Only 1 offset value can be entered. The entered offset value is valid for complete temperature range. This means if the setpoint temperature is changed the same offset value remains.

## 8. Cleaning

Make sure to remove the machine from the wall plug before cleaning and do not use abrasive cleaners. We suggest using a damp cloth.

## 9. Ordering Information

<b>Cat No.</b>	<b>Descriptions</b>
6.263 572	LLG-uni <i>INC</i> U 20 Cool Cooling Incubator, EU plug
6.263 573	LLG-uni <i>INC</i> U 20 Cool Cooling Incubator, UK plug

## 10. Warranty

This product is warranted to be free from defects in material and workmanship for a period of three (3) years from date of purchase.

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in this instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces or other causes not arising from defects in original material or workmanship. This warranty does not cover any incidental or consequential damages, commercial loss or any other damages from the use of this product.

The warranty does not cover damage to paint or finish and defects or damages caused by physical and chemical abuse or normal wear and tear. The warranty is invalidated by any nonfactory modification, which will immediately terminate all liabilities on us for the products or damages caused by its use. The buyer and its customer shall be responsible for the product or use of products as well as any supervision required for safety. If requested the products must be returned to the distributor in well packed and insured manner and all shipping charges must be paid.

Information on the disposal of electrical and electronic devices in the European Community:

The disposal of electrical devices is regulated within the European Community by national regulations based on the European Directive 2012/19/EU on waste electrical and electronic equipment (WEEE). According to these regulations, any devices supplied after 13.06.2005 in the business to business sphere, to which this product is assigned, may no longer be disposed of in municipal or domestic waste. They are marked with the following symbol to indicate this.

As disposal regulations within the EU may vary from country to country, please contact your supplier if necessary.

### PRODUCT DISPOSAL

In case the product is to be disposed of, the relevant legal regulations are to be followed.

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**Lab Logistics Group GmbH**