Operation Manual

uniSHAKER 2 3D Tumbling Shaker

Introduction:

LLG-uniSHAKER 2 is a tumbling (nutating) mixer with the angle and speed set for gentle, but thorough, no/low foam tube mixing and other nutating mixer applications. The Mixer comes with a platform designed to fit inside the LLG-uni/NCU 20 mini incubator.

I. Specifications:

Speed: Fixed 20 rpm
Tilt angle: +/- 20 ° (from the horizontal)
Motor: Long life AC gear motor
Platforms: 15.25 x 20.25 cm
Mat: Flat & Dimpled, autoclavable
Load Capacity: 1.6kg centered and balanced
Dimensions: 20.3 x 15.25 x 20.3 cm
Weight: 2 kg
Electrical: 230 VAC, 50 Hz, 0.2 A
Operating environment: 4 °C to 45 °C
Warranty: 3 years

II. Product Set-Up:

To install the LLG-uniSHAKER 2 please first remove the drip tray from the base of your LLG-uni/NCU 20 mini incubator (see figure A.). Then plug the power cord into the power inlet on the back wall of the incubator chamber (see figure B.) Install the shaker on the floor of the incubator (it should not be installed on any of the included shelves).

Once the shaker has been installed, center the base so that the platform will not contact the front door or the back fan cover when rotating.

Figure A.

Figure B.

III. Product Operation:

Load the platform from the center out, making sure the items placed on the platform form a balanced load around the center of the platform. This is important as larger, overhung loads may cause the rocker to stall or tip. Turn the unit on by activating the ON/OFF switch so that it is illuminated. The rotator will begin to rotate.

IV. Care and Maintenance:

No routine maintenance is required other than to keep the unit clean. Cleaning can be done with a damp cloth. Avoid the use of solvents as they may attack the rubber mat of plastic (ABS) housing.

Service and Contact:

In the event of a warranty or technical issue please contact **your local LLG-Labware distributor.**

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Lab Logistics Group GmbH